


# Setting up Element

## CONNECTING ELEMENT AND ELECTRODES

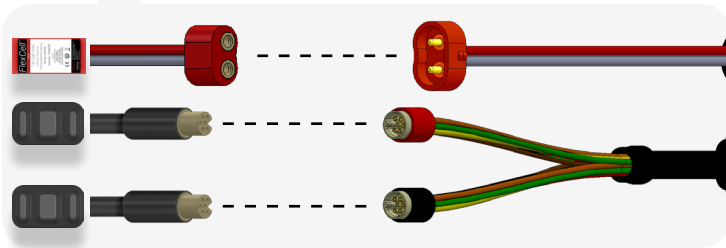


Consult the Instruction Manual for more details.

Standard 4-band coaxial plug shown.  
If your output connectors do not match, consult instruction manual.



And you're set!

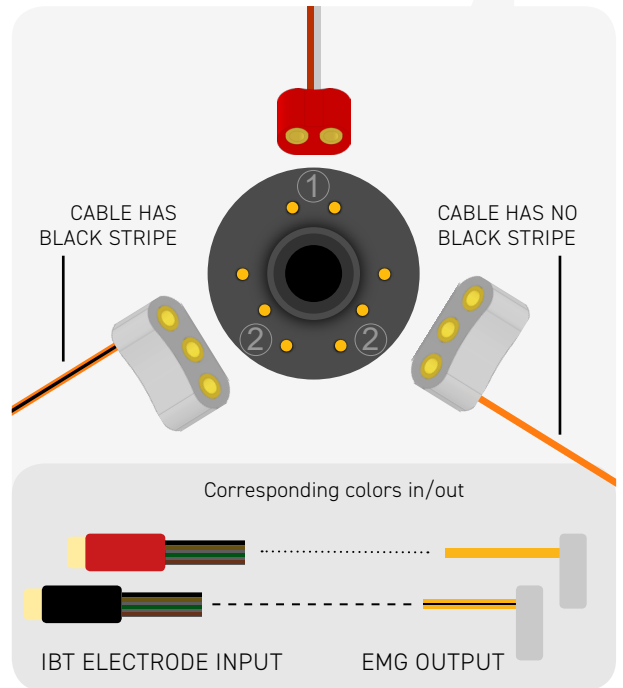

## CONNECTING ELEMENT TO PC OVER BLUETOOTH

Things You'll Need

PROVIDED IN BOX



Bluetooth Module

Just as a backup



Software Thumbdrive



Your PC

Minimum requirements:

Windows 10 (preferred) 500 MB Storage  
1.9 GHz Processor 1260 x 768 monitor  
4 GB Memory 1 USB Port



If your computer doesn't have it already

Installation instructions on next page

# Software Setup

- 1**

USB Paradox applies\*  
(See bottom)

Element Installer.msi  
Run me

Java Installer 32-bit.exe  
Java Installer 64-bit.exe

If needed  
(Choose 32 or 64-bit)

1.5  
If needed
- 2**

Start Menu

Type to search:  
"control panel"
- 3**

Hardware and Sound  
View devices and printers  
Add a device
- 4**

Add a device
- 5**

IBT Devices are always "IBTCore" and a 6-Digit number
- 6**

Launch Catalyst

Choose your Element device



And you're set! Enjoy!

## Troubleshooting Bluetooth

<b>ELEMENT CAN'T CONNECT</b>	<p>...to the Catalyst GUI after pairing</p> <p>...because I don't see its ID in Devices &amp; Printers</p>	<p>Power Element off then on. Element turns off its Bluetooth antenna if not connected for 10 minutes.</p>
<b>CATALYST: "BLUETOOTH NOT ACTIVE"</b>	<p>...even though I plugged the adapter in</p>	<p>Close then re-open Catalyst</p> <p>Unplug Bluetooth module and restart your computer, then plug in Bluetooth, wait 2-3 minutes, and open Catalyst</p>

If all else fails, call us at (443) 451-7175 and we'll help you out

\*USB Paradox (noun):  
The circumstance in which a USB plugs into the port only after attempting both sides